

For Research Use Only. NOT for Use in Diagnostic Procedures.

Extraction of *gamma*-Hydroxybutyric Acid (GHB) from Urine Using ISOLUTE® SLE+ Prior to GC/MS Analysis

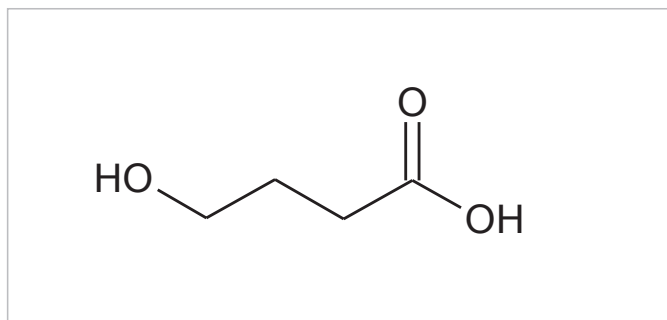


Figure 1. Structure of *gamma*-Hydroxybutyric acid (GHB)

Introduction

This application note describes the extraction of GHB from urine using supported liquid extraction and subsequent analysis by GC/MS.

ISOLUTE® SLE+ Supported Liquid Extraction plates and columns offer an efficient alternative to traditional liquid-liquid extraction (LLE) for bioanalytical sample preparation, providing high analyte recoveries, no emulsion formation, and significantly reduced sample preparation.

Analytes

Gamma-Hydroxybutyric acid (GHB) & GHB-D6

Sample Preparation Procedure

- Sample Pre-treatment:** Dilute pre-treated urine (0.2 mL) with 0.2% formic acid (aq) (0.2 mL). Spike GHB-D6 internal standard and vortex mix thoroughly.
- Format:** **ISOLUTE® SLE+ 200 µL Fixed Well Plate, part number 820-0200-P01**
- Sample Loading:** Load the pre-treated urine (200 µL total volume) onto each well and apply a pulse of vacuum or positive pressure (3–5 seconds) to initiate flow. Allow the sample to absorb for 5 minutes.
- Analyte Extraction:** Apply ethyl acetate (1 mL) and allow to flow under gravity for 5 minutes. Apply vacuum or positive pressure to pull through any remaining extraction solvent (5–10 seconds).
- Format:** **ISOLUTE® SLE+ 400 µL Sample Capacity columns, part number 820-0055-B**
- Sample Loading:** Load the pre-treated urine (200 µL total volume*) onto the column and apply a pulse of vacuum or positive pressure (3–5 seconds) to initiate flow. Allow the sample to absorb for 5 minutes.
- *Note:** Column is underloaded
- Analyte Extraction:** Apply ethyl acetate (1 mL) and allow to flow under gravity for 5 minutes. Apply a further aliquot of ethyl acetate (1 mL) and allow to flow for another 5 minutes under gravity. Apply vacuum or positive pressure to pull through any remaining extraction solvent (5–10 seconds).
- Post Elution, Reconstitution and Derivatisation:** Dry the extract in a stream of air or nitrogen at ambient temperature using a SPE Dry (20 to 40 L/min) or TurboVap (1.0 bar) for 30 mins.
- Upon dryness, reconstitute with ethyl acetate (50 µL) and BSTFA/1% TMCS (50 µL) and vortex for 20 seconds. Transfer to a high recovery glass vial.

GC Conditions

Instrument:	Agilent 7890A with QuickSwap
Column:	Phenomenex Zebron ZB-35, 30 m x 0.25 mm ID x 0.25 µm
Carrier	Helium 1.2 mL/min (constant flow)
Inlet:	250 °C, Split (ratio 10:1), 12 mL/min, septum purge flow: 3 mL/min
Injection:	1 µL
Wash Solvent:	Ethyl acetate
Oven:	Initial temperature 70 °C Ramp 10 °C/min to 140 °C Ramp 120 °C/min to 350 °C
Post Run:	Back-flush for 1.6 minutes (2 void volumes)
Transfer Line:	280 °C

MS Conditions

Instrument:	Agilent 5975C
Source:	230 °C
Quadrupole:	150 °C
MSD mode:	SIM

SIM Parameters

Table 1. Ions acquired in the selected Ion Monitoring (SIM) mode

SIM Group	Analyte	Quantifier Ion	Quantifier Ion
1	GHB-D6	239	240
1	GHB	233	117

Results

The optimized protocols for ISOLUTE SLE+ plates and columns demonstrated reproducible recoveries between three unique donors, as shown in Figure 2. Percentage recovery was 52–69%. RSDs for GHB were below 9% on the 200 µL capacity plate format and below 4% on 400 µL capacity column format. **Figure 3.** shows linearity data for plate and column formats ($r^2 > 0.999$ for both formats). This application is unusual because of the relatively high concentration cut-off requirement (10 µg/mL) for illicit GHB. The low matrix volume used and relatively low recoveries do not adversely affect the required assay performance and sufficient sensitivity was achieved.

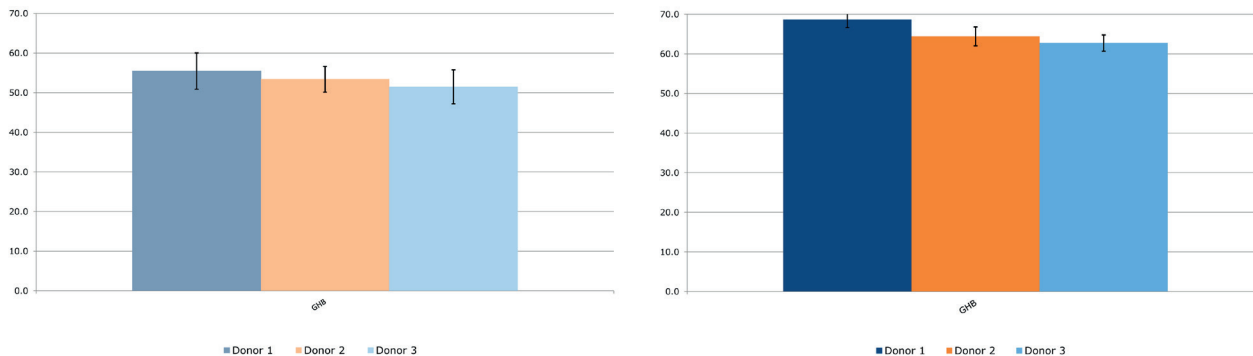


Figure 2. Typical analyte % extraction recoveries (n=7) using the ISOLUTE® SLE+ protocol using 200 µL capacity plate format (left) and 400 µL capacity column format (right)

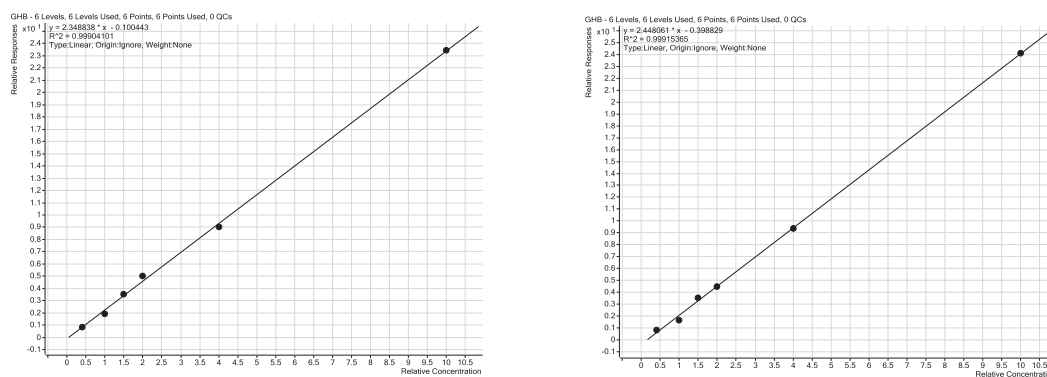


Figure 3. Calibration curves for extracted levels of spiked urine, using ISOLUTE® SLE+ protocol on 200 µL capacity plate format (left) and 400 µL capacity column format (right). Concentrations are 2.5, 5, 7.5, 10, 20 and 50 µg/mL. r^2 values are greater than 0.999.

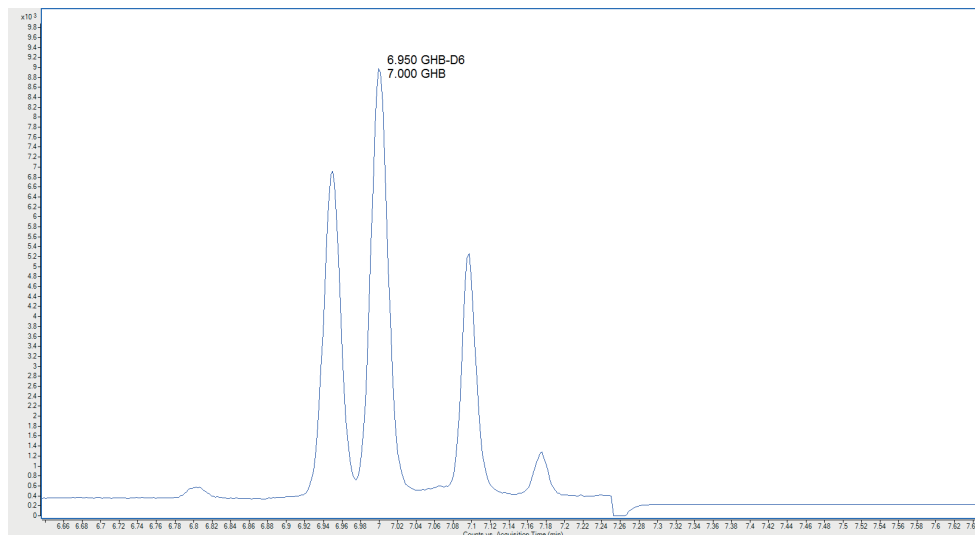


Figure 4. GC/MS chromatography for urine (Cal 1) spiked at 2.5 ng/mL. Compared to other more traditional solid phase extraction procedures, this ISOLUTE® SLE+ method results in a baseline with reduced noise and thus offers greater confidence when qualifying GHB.

Additional Information

0.2% formic acid (aq) is prepared by adding 100 µL concentrated formic acid (commercially available 98%) to 49.9 mL HPLC grade water

Ordering Information

Part Number	Description	Quantity
820-0200-P01	ISOLUTE® SLE+ 200 µL Supported Liquid Extraction Plate	1
820-0055-B	ISOLUTE® SLE+ 400 µL Sample Volume Columns	50
PPM-48	Biotage® PRESSURE+ 48 Positive Pressure Manifold	1
SD-9600-DHS-EU	Biotage® SPE Dry Sample Concentrator System 220/240 V	1
SD-9600-DHS-NA	Biotage® SPE Dry Sample Concentrator System 100/120 V	1
C103198	TurboVap® LV, 100/120V	1
C103199	TurboVap® LV, 220/240V	1

For the latest application notes visit www.biotage.com

EUROPE

Main Office: +46 18 565900
 Toll Free: +800 18 565710
 Fax: +46 18 591922
 Order Tel: +46 18 565710
 Order Fax: +46 18 565705
order@biotage.com
 Support Tel: +46 18 56 59 11
 Support Fax: +46 18 56 57 11
eu-1-pointsupport@biotage.com

NORTH & LATIN AMERICA

Main Office: +1 704 654 4900
 Toll Free: +1 800 446 4752
 Fax: +1 704 654 4917
 Order Tel: +1 704 654 4900
 Order Fax: +1 434 296 8217
ordermailbox@biotage.com
 Support Tel: +1 800 446 4752
 Outside US: +1 704 654 4900
us-1-pointsupport@biotage.com

JAPAN

Tel: +81 3 5627 3123
 Fax: +81 3 5627 3121
jp_order@biotage.com
jp-1-pointsupport@biotage.com

CHINA

Tel: +86 21 2898 6655
 Fax: +86 21 2898 6153
cn_order@biotage.com
cn-1-pointsupport@biotage.com

To locate a distributor,
 please visit our website at
www.biotage.com

Part Number: AN828.V.1

© 2014 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. Product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe. **FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.**

For more information visit www.biotage.com.