

For Research Use Only. NOT for Use in Diagnostic Procedures.

Extraction of Phencyclidine (PCP) from Urine Using ISOLUTE® SLE+ Prior to GC/MS Analysis

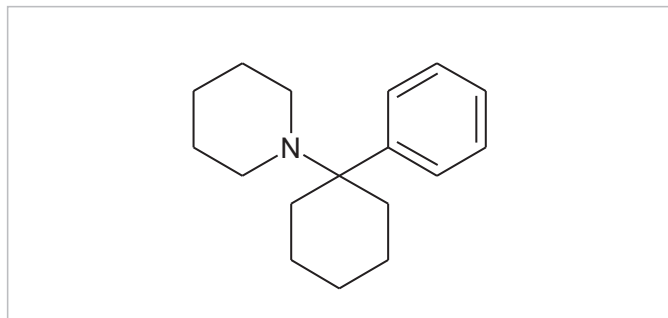


Figure 1. PCP structure

Introduction

This application note describes the extraction of PCP from urine using supported liquid extraction and subsequent analysis by GC/MS.

ISOLUTE® SLE+ Supported Liquid Extraction plates and columns offer an efficient alternative to traditional liquid-liquid extraction (LLE) for bioanalytical sample preparation, providing high analyte recoveries, no emulsion formation, and significantly reduced sample preparation.

Analytes

Phencyclidine (PCP) & PCP-D5

Sample Preparation Procedure

- Sample Pre-treatment:** Dilute pre-treated urine (1 mL) with 0.5% ammonium hydroxide (aq) (1 mL). Spike PCP-D5 internal standard and vortex mix thoroughly.
- Format:** **ISOLUTE® SLE+ 1 mL Sample Volume Columns, part number 820-0140-C (also available in tabless format, part number 820-0140-CG)**
- Sample Loading:** Load pre-treated urine (1 mL) onto the column and apply a pulse of vacuum or positive pressure (3–5 seconds) to initiate flow. Allow the sample to absorb for 5 minutes.
- Analyte Extraction:** Apply 1-chlorobutane (2.5 mL) and allow to flow under gravity for 5 minutes. Apply a further aliquot of 1-chlorobutane (2.5 mL) and allow to flow for another 5 minutes under gravity. Apply vacuum or positive pressure to pull through any remaining extraction solvent (5–10 seconds).
- Post Elution and Reconstitution:** To each sample, add 0.2M methanolic HCl (100 µL).
Dry the extract in a stream of air or nitrogen at ambient temperature using a SPE Dry (20 to 40 L/min) or TurboVap (1.0 bar) for 45 mins. Upon dryness, reconstitute with ethyl acetate (200 µL) and transfer to a high recovery glass vial.

GC Conditions

Instrument:	Agilent 7890A with QuickSwap
Column:	Agilent J&W DB-5 30 m x 0.25 mm ID x 0.25 µm
Carrier	Helium 1.2 mL/min
Inlet:	175 °C, Splitless, purge on at 2.0 minute, 50 mL/min
Injection:	1 µL
Wash Solvents:	1: Acetone 2: Ethyl acetate
Oven:	Initial temperature 60 °C, hold for 2 mins Ramp 80 °C/min to 200 °C, hold for 6.25 mins Ramp 100 °C/min to 280 °C
Post Run:	325 °C, Back-flush for 2.4 minutes (3 void volumes)
Transfer Line:	280 °C

MS Conditions

Instrument:	Agilent 5975C
Source:	230 °C
Quadrupole:	150 °C
MSD mode:	SIM

SIM Parameters

Table 1. Ions acquired in the selected Ion Monitoring (SIM) mode

SIM Group	Analyte	Quantifier Ion	Quantifier Ion 1	Quantifier Ion 2
1	PCP-D5	205	96	
1	PCP	200	91	130

Results

The optimized protocol demonstrated recoveries of greater than 88% from three unique urine donors, with corresponding RSDs of less than 8% (n=7), illustrated in **Figure 2**. **Figure 3**. shows the calibration line formed from the extraction of spiked urine at levels of 10 ng/mL up to 1000 ng/mL.

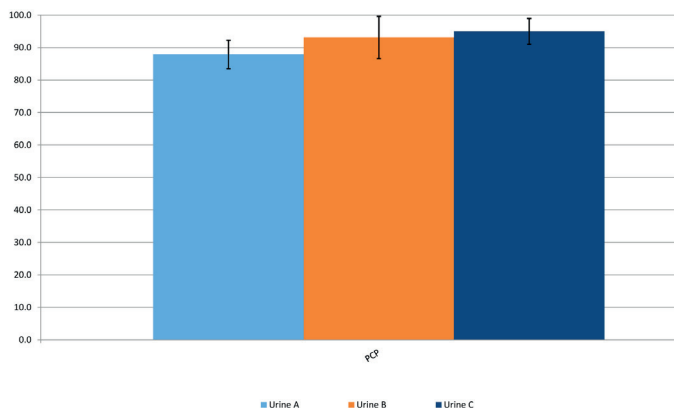


Figure 2. Typical analyte % extraction recoveries (n=7) using the ISOLUTE® SLE+ protocol on 1 mL C column.

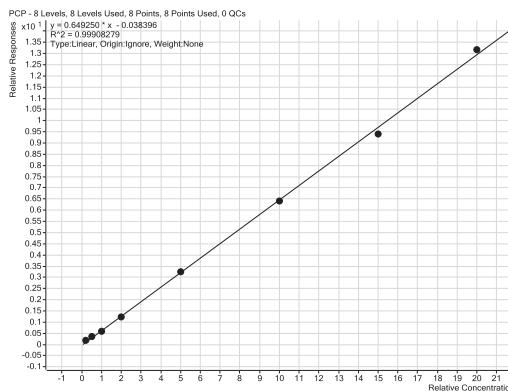


Figure 3. Calibration line for extracted levels of spiked urine, using ISOLUTE® SLE+ protocol described. The coefficient of determination was determined to be greater than 0.999 across concentration values 10, 25, 50, 100, 250, 500, 750 and 1000 ng/mL. The lower limit of quantitation (LLOQ) was determined to be 10 ng/mL.

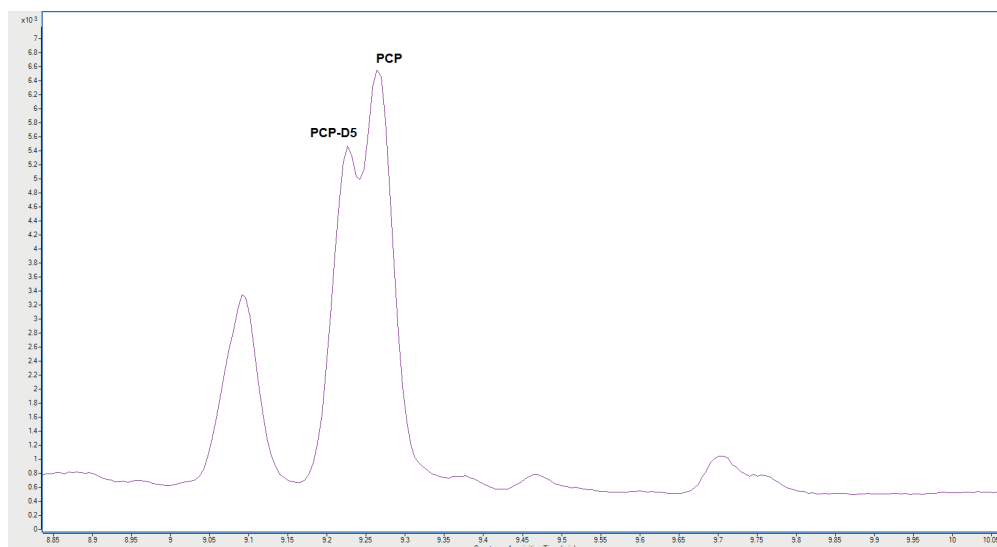


Figure 4. GC/MS chromatography for urine spiked with both PCP and PCP-D5 at 100 ng/mL. TIC acquisition of m/z 205, 96, 200, 91, 130

Additional Information

0.5% ammonium hydroxide (aq) is prepared by adding 0.5 mL concentrated ammonium hydroxide (commercially available 28%–30% NH₃ in H₂O) to 99.5 mL HPLC grade water.

0.2M methanolic HCl is prepared by adding 200 µL concentrated HCl (commercially available 37%) to 11.8 mL HPLC grade methanol.

Ordering Information

Part Number	Description	Quantity
820-0140-C	ISOLUTE SLE+ 1mL Supported Liquid Extraction Column	30
820-0140-CG	ISOLUTE SLE+ 1mL Supported Liquid Extraction Column (Tablets)	30
PPM-48	Biotage® PRESSURE+ 48 Positive Pressure Manifold	1
SD-9600-DHS-EU	Biotage® SPE Dry Sample Concentrator System 220/240 V	1
SD-9600-DHS-NA	Biotage® SPE Dry Sample Concentrator System 100/120 V	1
C103198	TurboVap® LV, 100/120V	1
C103199	TurboVap® LV, 220/240V	1

For the latest application notes visit www.biotage.com

EUROPE

Main Office: +46 18 565900
 Toll Free: +800 18 565710
 Fax: +46 18 591922
 Order Tel: +46 18 565710
 Order Fax: +46 18 565705
order@biotage.com
 Support Tel: +46 18 56 59 11
 Support Fax: +46 18 56 57 11
eu-1-pointsupport@biotage.com

NORTH & LATIN AMERICA

Main Office: +1 704 654 4900
 Toll Free: +1 800 446 4752
 Fax: +1 704 654 4917
 Order Tel: +1 704 654 4900
 Order Fax: +1 434 296 8217
ordermailbox@biotage.com
 Support Tel: +1 800 446 4752
 Outside US: +1 704 654 4900
us-1-pointsupport@biotage.com

JAPAN

Tel: +81 3 5627 3123
 Fax: +81 3 5627 3121
jp_order@biotage.com
jp-1-pointsupport@biotage.com

CHINA

Tel: +86 21 2898 6655
 Fax: +86 21 2898 6153
cn_order@biotage.com
cn-1-pointsupport@biotage.com

To locate a distributor,
 please visit our website at
www.biotage.com

Part Number: AN829.V.1

© 2014 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. Product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe. **FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.**

For more information visit www.biotage.com.