


Biotage® PS-TsNHNH₂

Electrophile Scavenger

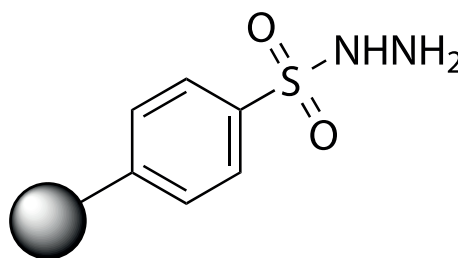

Key Facts



Shelf Life

Capacity
(mmol/g)

BSE/TSE



Scalable

Particle Size
(μm)Thermally &
Mechanically
StableGood
Laboratory
PracticeBulk Density
(g/L)

Specifications

| | |
|-------------------------------|---|
| Chemical Name: | Polystyrene sulfonyl hydrazide |
| Resin Type: | 1% Cross-linked poly(styrene-co-divinylbenzene) |
| Application: | Scavenging aldehydes and ketones |
| Scavenging Conditions: | 3 equivalents relative to carbonyl, 1–3 h, 20 °C, DCM. Ketones and hindered aldehydes are accelerated by the addition of HOAc (~10%) and/or heat. |
| Compatible Solvents: | DCM (7 mL/g), DCE (7 mL/g), THF (6.5 mL/g), DMF (7.2 mL/g) |
| Storage: | Cool, dry location |

PS-TsNHNH₂ is a resin-bound equivalent of p-toluenesulfonyl hydrazide and is an excellent scavenger of aldehydes and ketones. Comparison with other polymeric benzyl hydrazide showed PS-TsNHNH₂ was a superior scavenger for carbonyls and much more stable to storage (the benzyl hydrazide resin decomposed on storage).


Removal of excess carbonyls from solution generally requires a three-fold excess of PS-TsNHNH₂. Addition of a catalytic

amount of HOAc (5–10%) may be required for ketones and hindered aldehydes. HOAc is also required for sequestering aldehydes in DMF. Complete removal of common aldehydes occurs in 0.5 to 3 h and removal of a ketone takes from 2 to 16 h. Elevated temperatures were required for hindered ketones (e.g. 2,6-dimethylcyclohexanone). Upon completion of the scavenging, the resin is rinsed with a suitable solvent (i.e. those which swell polystyrene), and the product is isolated by concentration. PS-TsNHNH₂ was successfully used to work up the synthesis of alcohols by addition of a Grignard reagent to aldehydes.

PS-TsNHNH₂ is also potentially useful as a polymer-bound reagent. Bound sulfonyl hydrazones, formed by condensation with carbonyl compounds, can be utilized in further synthetic transformations. The high accessibility of tosyl hydrazide functional groups in PS-TsNHNH₂ afford high synthetic fidelity relative to reported systems.¹

Representative Procedure

PS-Tosyl Hydrazide (0.6 g) was added to a mixture of 3-methoxybenzaldehyde (20 μL, 1.5 mmol) and naphthalene (15 mg; internal standard) in THF (2 mL). This mixture was stirred for 30 min, then filtered and washed with MeOH (8 mL). The filtrate was concentrated and analyzed by RP-HPLC.

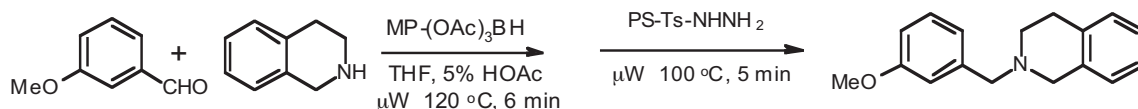

Scheme 2. One pot microwave-assisted reductive amination reaction followed by aldehyde scavenging.

Reductive Amination

This scavenger can be applied to the rapid clean up of tertiary amines prepared through reductive amination using MP-Triacetoxyborohydride and microwave irradiation (Scheme 1).

To a solution of the amine (0.5 mmol) and 3-methoxybenzaldehyde (0.7 mmol) in a solution of THF and 5 % Acetic acid (2 mL) was added MP-Triacetoxyborohydride (523 mg, 2.39 mmol/g, 1.25 mmol). The vial was capped and heated to 120 °C for 6 min, then PS-TsNHNH₂ (0.6 g) added. The vial mixture was heated to 100 °C for 5 min, after which the reaction mixture was filtered. The solid was washed with a 1:1 solution of THF/MeOH (10 mL) and the filtrate was concentrated to afford product.

Ordering Information

| Part Number | Quantity |
|-------------|----------|
| 800497 | 3 g |
| 800270 | 10 g |
| 800271 | 5 g |
| 800272 | 100 g |
| 800317 | 1000 g |

EUROPE

Main Office: +46 18 565900
 Fax: +46 18 591922
 Order Tel: +46 18 565710
 Order Fax: +46 18 565705
 order@biotage.com
 Support Tel: +46 18 56 59 11
 Support Fax: +46 18 56 57 11
 eu-1-pointsupport@biotage.com

NORTH & LATIN AMERICA

Main Office: +1 704 654 4900
 Toll Free: +1 800 446 4752
 Fax: +1 704 654 4917
 Order Tel: +1 800 446 4752
 Order Fax: +1 704 654 4917
 ordermailbox@biotage.com
 Support Tel: +1 800 446 4752
 us-1-pointsupport@biotage.com

JAPAN

Tel: +81 3 5627 3123
 Fax: +81 3 5627 3121
 jp_order@biotage.com
 jp-1-pointsupport@biotage.com

CHINA

Tel: +86 21 68162810
 Fax: +86 21 68162829
 cn_order@biotage.com
 cn-1-pointsupport@biotage.com

KOREA

Tel: +82 31 706 8500
 Fax: +82 31 706 8510
 korea_info@biotage.com
 kr-1-pointsupport@biotage.com

INDIA

Tel: +91 11 45653772
 india@biotage.com

Distributors in other regions are listed on www.biotage.com

Part Number: PPS403.V.2

© 2023 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. A list of all trademarks owned by Biotage AB is available at www.biotage.com/legal. Other product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe.